Crime-lite® ML2

Bench top magnifier with multi-wavelength illumination for general search and fluorescence examinations

- Optimised for the examination of large items in the laboratory
- Choice of UV, Violet, Blue, Blue/Green, Green, Orange, Red & IR LED modules
- For the fast and effective visualisation and magnification of:

Blood Fingerprints Body fluids Gunshot residues Accelerants Fibres etc.

foster + freeman

ASORATORY C

Crime-lite[®] ML2

BENCH TOP MAGNIFIER WITH MULTI-WAVELENGTH ILLUMINATION

The new **Crime-lite ML2** combines the versatility of multi-wavelength, high intensity LED illumination with wide area bi-ocular magnification to provide an effective bench top search and examination tool in the forensic laboratory.

The **Crime-lite ML2** features 8 high intensity white LEDs for general search and a choice of up to 4 additional narrowband LEDs for the examination of fluorescent evidence including:

UV, Violet, Blue, Blue/Green, Green, Orange, Red & IR

Designed for the examination and 'marking up' of large items of evidence in the laboratory, the **Crime-lite ML2** head is mounted on an extendable articulated arm with a reach of over 1 metre and includes a 3.5 dioptre lens that provides a nominal magnification of 1.8x.



The optional 5MP UV-Vis-IR camera is rotated into position over the magnifier lens and can be retro-fitted to existing Crime-lite ML light sources as a quick and easy image capture device.



The range of examination wavelengths has now been extended with the addition of a new Infrared module.

Using an UV-Vis IR sensitive camera (pictured left) the new module is capable of detecting blood, trace evidence and fibres on difficult backgrounds through the visualisation of infrared images.

Operation is via a PC or laptop loaded with image capture software.

Crime-lite[®] ML2 Applications

The Crime-lite ML2 provides forensic examiners with a fast and efficient method of screening articles for the presence of fingerprints, trace evidence or other biological material. Laboratory productivity is greatly increased by quickly 'marking up' areas of interest and eliminating articles devoid of evidence.

Use the look-up tabels below to determine the wavelengths required for your chosen applications...


	W	UV	V	В	BG	G	0	R	IR
Blood	\checkmark		\checkmark	\checkmark					\checkmark
Body Fluids		\checkmark	\checkmark	\checkmark	\checkmark				
Bites & Bruising		\checkmark	\checkmark						
Bone & Tooth				\checkmark	\checkmark	\checkmark			
Hair			\checkmark	\checkmark					

Biological & Trace Evidence

Fibres		\checkmark	\checkmark	\checkmark	\checkmark	\checkmark	\checkmark		\checkmark
GunShot Residues					\checkmark	\checkmark			\checkmark
Drug Residues		\checkmark	\checkmark						
Accelerants				\checkmark	\checkmark	\checkmark			
Shoeprints	\checkmark								
General Debris	\checkmark		\checkmark	\checkmark	\checkmark	\checkmark	\checkmark	\checkmark	

Fingerprints


	UV	V	В	BG	G	0	R	IR
Acid Yellow			\checkmark					
Ardrox			\checkmark					
NEW ASP Powder							\checkmark	\checkmark
Basic Yellow		\checkmark	\checkmark					
DFO				\checkmark	\checkmark	\checkmark		
Greenwop				\checkmark				
Magnetic Red			\checkmark					
Nanoparticles	\checkmark	\checkmark	\checkmark	\checkmark	\checkmark			
Ninhydrin					\checkmark	\checkmark	\checkmark	
PolyCyano UV	\checkmark							
Redwop			√	\checkmark				
Rhodamine 6G				\checkmark				

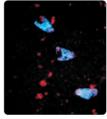

Human teeth



Fingerprints in blood

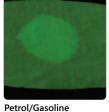












Redwop

Hair

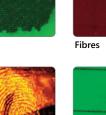







Ninhydrin





PolyCyano UV


Searching Beyond the Visible


The addition of a new Infrared LED module and UV-Vis-IR camera has further increased the functionality of the Crime-lite ML2 by allowing investigators to conduct examinations beyond the limits of the visible spectrum.

Previously hard to detect evidence, such as blood or GSR on dark or densely patterned fabrics, becomes clearly visible under IR illumination when viewed in the infrared waveband.




Blood on black wool

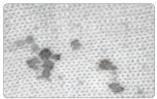

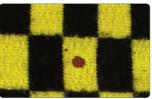

Blood on multi-coloured cotton

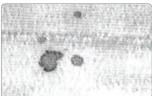

Crime-lite ML2 Specifications

The versatility of the design allows up to 4 narrow band LED modules to be fitted, in addition to a white light LED module.

CRIME-LITE ML2 Standard Features

White LED Module

ML2 Body

- 8 x High efficiency White LEDs
- 400-700nm (nominal)
- Adjustable output intensity
- Homogenous light beam •
- Large 3.5 dioptre magnifying lens Easy clean touch panel controls •
- Dust cover .
- Power input 100-240V, 80W
- Built-in thermal protection
- Heavy duty articulated arm
- Reach 132cm
- Bench or wall mounted
- 3-axis adjustable head mount • •
- Internal cables (no trailing wires)

Laboratory Arm

- Heavy duty articulated arm .
- Reach 132cm
- Bench or wall mounted .
- 3-axis adjustable head mount Internal cables (no trailing wires)

Narrow band LED Options


	UV module	Violet module	Blue module	Blue/Green module	Green module	Orange module	Red module	IR module
Number of LEDs	16 X	4 X	16 X	16 X	16 X	16 X	16 X	16 X
Peak bandwidth	365nm	405nm	445nm	475nm	520nm	590nm	640nm	860nm
10% bandwidth	350-380nm	395-425nm	420-470nm	445-510nm	480-560nm	570-610nm	600-660nm	800-900nm
Viewing filter	GG420	GG455nm	GG495nm	OG550nm	OG590nm	RG645nm	IR/Vis	IR/Vis

Custom Configurations

In addition to White LEDs, the Crime-lite ML2 can be fitted with up to four narrowband LED modules. Use the table on the right to select a valid combination.*


Narrowband LED modules are supplied together with the corresponding fluorescence viewing filters.

IR and Red LED modules require the UV-Vis-IR sensitive camera and connection to a PC running the Crime-lite Cam software (see next page for details)

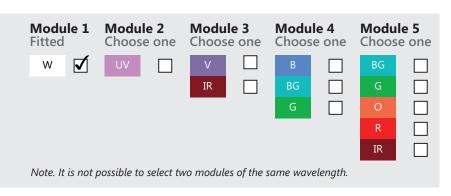

Common Configurations

The look-up table below includes details of the most commonly requested combinations.

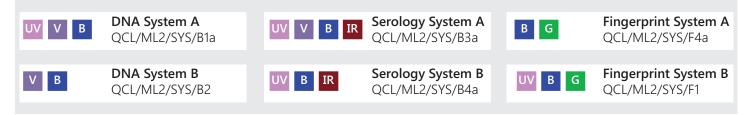

* Illumination module options must be specified and fitted during the equipments manufacture and cannot be retro-fitted to an existing unit.

If you have any questions regarding module selection or the Crime-lite ML2 please contact sales@fosterfreeman.com for further guidance.

Crime-lite[®] ML2 Accessories

QCL/ML2/CAM

QCL/ML2/PENT

QCL/ML2/PENT/LT

UV-VIS-IR Camera

Attaches to Crime-lite ML2 for and enables image capture

- 5MP CMOS auto-focus camera .
- Essential for the visualisation of IR images •
- 8mm f/1.8 lens .
- . Sensitive from 350nm to 1100nm
- Swings into position over viewing lens .
- . USB 2.0 interface
- Wireless remote control .
- 4-position filter wheel with UV, Vis & IR filters .
- Image capture software •

Computer systems

For use with VIS/IR Camera module.

PENTIUM COMPUTER

- High specification PC .
- 22" Flat screen TFT monitor

LAPTOP COMPUTER

- High specification laptop PC
- 17" monitor

Contact Foster+Freeman for current specifications

Contrast enhancing filters

For use with White LEDs

BLUE	for blood search	QCL/ML/CONT/B
GREEN	for viewing exhibits treated with Ninhydrin	QCL/ML/CONT/G

Bandpass viewing filters

Tailored for the examination of specific evidence types

QCL/ML/VFBP415 415nm visualisation of blood spatter (centered at 415nm) QCL/ML/VFBP530 530nm visualisation of body fluids (centered at 530nm) 550nm visualisation of body fluids (centered at 550nm) QCL/ML/VFBP550

Laboratory Cart

For easy positioning and high mobility

 LABORATORY CART Wheeled cart with mounting plate for the articulated arm plus UPS storage space 	QCL/245
ADJUSTABLE LAPTOP MOUNTFor use with Laboratory Cart	QCL/247
UPS (not shown) Run time 4 hours 	QCL/246

Consumables

PARTICULATE AIR FILTER • Replacements for the ML2, Pack of 50 QCL/ML/FILTPACK









* All specified LED wavelengths are nominal. Foster + Freeman reserve the right to alter the specification of the Crime-lite ML2 and its accessories without prior notice.



The Crime-lite range

Crime-lite[®] XL

The worlds most powerful handheld LED forensic light source, the Crime-lite XL features 96 surface mount LEDs providing up to 115 Watts of radiant power capable of revealing the minutest traces of evidence.

Crime-lite[®] 825

Where serious crime demands intensive crime scene investigation, the 16 LED Crime-lite 82S offers high levels of illumination.

Crime-lite[®] 82L

16 high intensity white LEDs mounted behind a cylindrical lens give a wide, oblique light beam that is ideal for detecting surface debris and shoe prints in dust.

Crime-lite[®] 425

The Crime-lite 42S provides three illumination outputs in a single light source.

Crime-lite[®] 2

Effective illumination of evidence from a single high intensity LED.

WHY CHOOSE CRIME-LITE®?

- High quality of design and construction
- LEDs selected for peak performance and wavelength
- 100% light output throughout battery charge
- Worldwide product support
- · High power to weight ratios
- · Compact and energy efficient

	No. of LEDs	LED output po	No. of Wavele	Rechargeable	A/C Mains Op	Weight
Crime-lite 2	1	1.2W	8	Yes	Yes	490g
Crime-lite 82S	16	19W	8	Yes	Yes	560g
Crime-lite 82L	16	N/A	8	Yes	Yes	900g
Crime-lite 42S	8+8	9W	6	Yes	Yes	560g
Crime-lite XL	96	115W	4	Yes	Yes	1.75Kg

wer

Battery

ion

Head Office, UK Sales Office Vale Park | Evesham | WR11 1TD | United Kingdom

Tel: +44 (0)1386 768 050 | sales@fosterfreeman.com

USA Sales Office

46030 Manekin Plaza | Suite 170 | Sterling | VA 20166 | USA Tel: 888 445 5048 | usoffice@fosterfreeman.com

foster + freeman

fosterfreeman.com